

## Quick Guide Pipette Testing Procedure

- 1 Turn on ATMOS and select "Accuracy" on the main screen.
- 2 Select your pipette size.
- 3 Insert the pipette tip into the port.  
**NO LIQUID IN THE TIP!**
- 4 Press the button **Start** on the screen.
- 5 Press the piston down when the "Press and Hold" image appears on screen.  
The displacement should be done in about one second (**blue progress bar**).  
Don't move over the first stop of the piston.
- 6 Keep stable during the hold period (**red progress bar**).
- 7 Release the piston when "Release" appears on the screen (**green progress bar**). No need to remove the tip from the port until the next measurement.
- 8 Choose what to do with your measurement:
  - Skip: discard measurement, not stored in memory.
  - Next: start new measurement, current data is stored.
  - End: Stop series and show statistics (min. 3 measurements necessary).
- 9 After selecting "End", you see the average of your measurement series and the standard deviation. Select "New measurement" to start a new series or select another function.



Detailed manual:



[www.unipix-atmos.com/downloads](http://www.unipix-atmos.com/downloads)